

16 May 2004

How to Start cryo control window

Cryogenics control window is open on PC0 (near phone). If it is not, then start it as follows

1. Open Iexplorer and connect to <http://137.138.65.62/anibus.html>
2. Login as user [atlas](#) with psw [cern2](#)
3. Answer [Yes](#) to warning
4. Click [OK](#) (do not need password)
5. Click [Connect to CERN](#) and wait a few seconds when pictures apper
6. Click [Exit](#)
7. Click on [Main View](#)
8. Click on [HEC temperature](#)